


Multipette® / Repeater®

Caution! Immediate permanent sensor damage can occur.



- ▶ Never rotate the adapter when inserting!
- ▶ Never rotate the adapter once inserted!
- ▶ Never rotate any Combitip plus once inserted!
- ▶ For further information see manual.

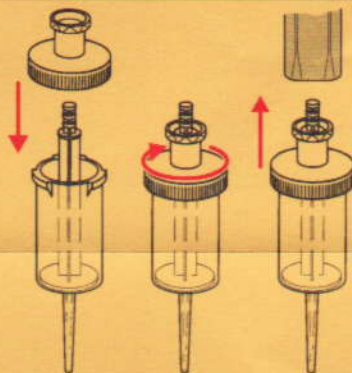

- ▶ Always connect the adapter to the Combitip plus before attaching to the Multipette / Repeater!

- ▶ Always hold the Multipette / Repeater by the housing.
- ▶ When using the 25 mL or 50 mL Combitips plus, ensure that the adapter is in perfect condition.
- ▶ The adapter is a wear part and has to be changed on a regular basis.

eppendorf

In touch with life

Your local distributor: www.eppendorf.com/worldwide

Eppendorf AG · 22331 Hamburg · Germany · Tel. +49 40 538 01-0
Fax +49 40 538 01-556 · E-Mail: eppendorf@eppendorf.com

Eppendorf North America, Inc. · One Cantiague Road, P.O. Box 1019
Westbury, N.Y. 11590-0207 USA · Tel. +1 516 334 7500
Toll free phone 800 645 3050 · Fax +1 516 334 7506
E-Mail: info@eppendorf.com

Application Support

Europe, International: Tel. +49 1803 666 789 · E-Mail: support@eppendorf.com
North America: Tel. 800 645 3050 ext. 2258 · E-Mail: support_NA@eppendorf.com
Asia, Pacific: Tel. +603 8023 2769 · E-Mail: support_Asia@eppendorf.com

eppendorf® is a registered trademark. 0012 573 634-00/0807